



Laboperator v3.4.1

Changelog

2021-07-06

---

## Changelog

---

### v3.4.1 - 2021-07-06

---

#### Changed

- [Dashboards] When adding Gauge and Gauge Control elements into a dashboard, the respective min/max value will now be automatically populated from the min/max value defined by the related device's channel
- [Workflows] Minor UX change in the Change Reason modal on which the field value will show up above the reason.

#### Fixed

- [Collections] Minor bug fixes and UX improvements
- [Workflows] Timer which are started in one substep can now trigger completion event in subsequent steps and substeps
- [Workflows] Performance improvement when creating workflow runs with a lot of fields

### v3.4.0 - 2021-06-22

---

#### Added

- [Audit Trails] Enable exports of the audit trails into csv/pdf/xlsx format. This feature depends on the report engine module which will be generally available soon. For on-premise installation, contact us for support if you want to have access before it is generally available.
- [Collections] Allow filtering collection resources by role. This can be useful when reviewing the roles and rights granted to members of the organization.
- [Workflows] Add new updateResource action which can be used to update related resources from within a running workflow. For now, this action is limited to updating the title of the running workflow.

#### Changed

- [Collections] Improve accessibility of various status indicators.
- [Audit Trails] Change format of date time picker to be more GxP compliance.

## Fixed

- [Notifications] Fix delivery of notifications via email
- [Users] Fix redirect after user sign in via SSO. Users will now correctly be redirected to the page they were trying to access before the authentication.
- [Webhooks] Fix display of headers when editing webhook subscriptions
- [Workflows] Fix potentially data lost when updating workflow tables in very quick succession under a slow network.

## v3.3.1 (Hotfix) - 2021-06-10

---

### Fixed

- [Collections] Fix page freezing on certain resource navigation.

## v3.3.0 - 2021-06-10

---

### Added

- [Collections] Organisation administrators will now be able to navigate to resources that other members have access to.
- [Collections] Provide ways for users to review their roles across multiple resources

### Fixed

- [Audit Trails] [Workflows] Fix invalid translations and missing validation messages.

## v3.2.0 - 2021-05-27

---

### Added

- [Audit Trail] Add support for filtering based on time range and increase the allowed page size

## Changed

- [Exports] Export to network storage is now deprecated and will be removed with the next major release. The corresponding feature flag within the organization setting is now frozen.
- [Workflows] Actions which are triggered in short succession during a workflow run will now deterministically be resolved in the sequence of when the action is triggered. Which should considerably improve the experience of working with complicated workflows

## Fixed

- [Devices] Fix broken file upload on data point for certain file types
- [Workflows] Fix support for reordering upcoming steps

## v3.1.1 - 2021-05-13

---

## Fixed

- [Audit Trail] Various bug fixes related to entry display
- [Dashboards] Fix display when some SI units is used by the device
- [Devices] Fix device setting page not showing when the device name contains Unicode characters
- [Workflows] Various bug fixes related to timer behaviors

## v3.1.0 - 2021-04-29

---

## Added

- [Audit Trail] Improve UI/UX for discovering and filtering related application events

## Fixed

- [Devices] Various bug fixes and performance improvements
- [Workflows] Various bug fixes and performance improvements

## v3.0.1 - 2021-04-15

---

## Fixed

- [Collections] Organisation admin can now correctly archive all resources within the organisation.
- [Workflows] In some cases, updates to fields which should have required change reason were accepted even when the change reason is missing. Such updates will be rejected as expected
- [Workflows] Scan pattern for each steps and substeps are now correctly detected as workflow executor progress to the steps
- [Workflows] Various other bug fixes and performance improvements

## v3.0.0 - 2021-04-01

### Added

- [Versions] UI/UX improvement on activities display

### Changed

- [Audit Trail] The Activities page is revamped with UI/UX improvements.

### Fixed

- [Connector][Dashboards] Improve rendering performance of some elements
- [Workflows] Fix various UI eligibility issue

### Removed

- [Job] Jobs and Tasks feature are removed
- [Deployment & Maintenance] Subdomain-based tenant strategy is removed in favour of route-based tenant strategy