



Laboperator v4.0.0

Changelog

2022-05-20

Changelog

v4.0.0 - 2022-05-20

Added

- [Devices] UR10 from Universal Robots: new device (driver-ur-dashboard v1.0.0).

Changed

- [Resources] Archival of resources are replaced with hiding of resources. The changes are mostly cosmetic with minimum expected impact on user's ability to manage their resources.
- [Devices] R-300, R-220 pro, SyncorePlus from Büchi: simplified driver parameters. An upgrade to this version requires manually updating the parameters. (driver-buchi-openinterface v2.0.0).

Removed

- [API] APIv1 of Laboperator was deprecated since release of Laboperator v2.17.0 and is finally removed
- [Administrations] Management of user permissions and access to resources via ACL is removed in favor of RBAC.
- [Audit Trail] Activities page is removed in favor of Audit Trail
- [Collections] With the removal of ACL, the ability to review ACL permissions and filter resources based on ACL roles is also removed
- [Connectors] Non-functional attributes are removed from the Connector edit form and info panel.
- [Exports] Support for exporting to remote Network Drive is removed.
- [Organizations] The explicit distinction between Administrator and Member is removed in favor of assigned roles and permissions via RBAC

Fixed

- [Dashboards] Fixed issue with invalid count of data points in Line Chart elements
- [Exports] Fixed issue with export with sufficiently large amount of data points
- [Workflows] Fixed issue with outdated snapshot observed in workflows which have behaviors written to sequentially update a field and complete substep/step within a single action.